

Ingram-Ward
Petrographic
Thin Section Cut-Off Saw
Models 135 and 137U

User's Manual


v. 1.88

Ward's Natural Science Establishment, Inc.

CONTENTS

Introduction	1
Unpacking Your Equipment	1
Setting Up	
Coolant System	3
Vacuum System	4
Electrical Hookup	6
General Operating Procedures	
Saw Blade Preparation	7
Making Thin Sections	7
Cut-Off Saw Operation	8
Maintenance	10
Troubleshooting	11
Keys to Better Thin Section Preparation	
Introduction/What is a Thin Section?	12
Specimen Preparation	13
Grinding the Chip	14
Microscope Slide Variations	16
Chip Mounting	16
Final Preparation	18
Customer Support	19
Support Equipment	20

Ingram-Ward Petrographic Thin Section Cut-Off Saw Models 135 and 137U

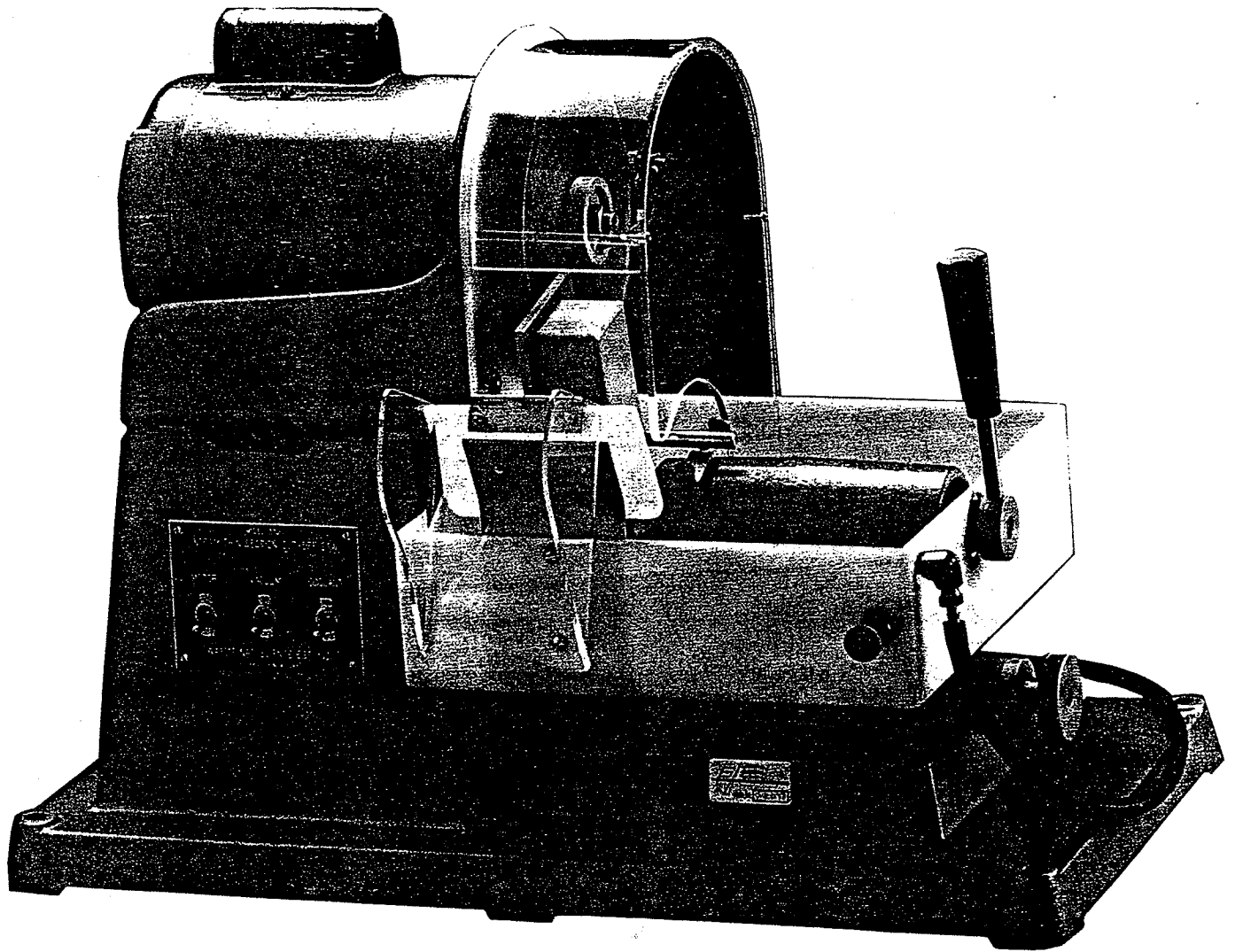

INTRODUCTION

This manual will familiarize you with the basic set-up and maintenance of the Ingram-Ward Petrographic Thin Section Cut-Off Saws, Models 135 and 137U, and present some recommended procedures and techniques for producing quality thin sections.

UNPACKING YOUR EQUIPMENT

Your Ingram-Ward Thin Section equipment has been packed with the utmost care for damage-free shipping. If, however, damage has occurred during shipment, please notify Ward's Customer Service Department and retain all crates and packing materials for inspection.

To unpack your equipment:

- (1) Remove the shipping crate top by taking out all perimeter fasteners.
- (2) Remove the 2" x 4" stud above the machine by removing the fasteners on the sides of the shipping crate. Remove any items, such as dressing sticks, rubber pads, etc., which may be attached to the stud.
- (3) Remove nuts and washers, found at the four corners of the machine base, which secure the unit to the bottom of the shipping crate.
- (4) Check to see that the knots on the lifting ropes are tight. Using these ropes, slowly lift the machine from the shipping crate, being sure to keep the machine level. This usually requires two people. The motor end is much heavier than the opposite end, so go slowly.
- (5) Place the machine on a sturdy table of proper working height (36 inches is recommended). An optional mobile lab table is also available from Ward's for more portable use. See your Ward's Geology/Earth Science Catalog for details.
- (6) Remove the lifting ropes from under the machine and place a rubber pad (provided) under each corner. CAUTION: DO NOT fasten the machines to the table. Doing so may distort the base and casting and cause damage to the equipment.

With your equipment package, you should find the following items:

- 1 Ingram-Ward Petrographic Thin Section Cut-Off Saw
(Model 135 or 137U)
- 4 Rubber Pads
- 1 Dressing stick for saw blade
- 3 Types of tubing for vacuum pump, coolant pump, and drain

Should any parts be damaged or missing, please contact Ward's Customer Service immediately for replacement.

NOTE:

Operation of this equipment requires the use of a coolant system and vacuum pump. These may be supplied by the user, or they are available separately from Ward's. Consult your Ward's catalog for details.

SETTING UP

Coolant System


Ingram-Ward Thin Section Equipment can be used with a self-contained coolant pump system (available separately) or the equipment may be connected directly to a water line.

- (1) **Coolant Pump System (Ward's)** -- Remove the plastic cover from the coolant pump reservoir. Place the recirculating pump in the bottom of the reservoir. Connect the clear plastic tubing (1/2" diameter) directly to the pump nozzle or to the connecting tube in the rubber stopper and through the plastic top. Connect the other end to the water supply intake at the rear of the thin section machine. Connect the larger plastic drain tubing (3/4" diameter) to the fitting at the back of the water catch pan (white plastic pan on machine) and route the other end to the filtered intake on the reservoir tank. Fill the coolant reservoir at least three quarters full with water or other desired coolant. (CAUTION: kerosene, toluene, or other solvent-based coolants may damage reservoir and water catch pan. Avoid their use.) Check and maintain coolant level periodically. Drain and clean sludge from the coolant pump reservoir as needed. This is important, since accumulated sludge can be recirculated by the pump, causing scratching of the microscope slide, the thin section, and the plastic coolant hood. Check to make sure all water line connections are secure.
- (2) **Direct Water Line** -- If the Ingram-Ward equipment is to be connected to a direct water line, be sure to install a shut-off valve in the water line before the nozzle connection on the machine. This shut-off valve is extremely important, as the plastic tubing will not hold the static pressure of most water systems. DO NOT open the shut-off valve completely. This valve should be used to adjust the initial water flow. Since this water will not be recirculated, the coolant return tubing should be connected to the fitting at the base of the plastic catch pan and directed to an existing drain. Any drain used should have a trap that can be accessed for the removal of sediment buildup. Check the drain trap periodically for sludge accumulation and clean as needed to avoid drain stoppage.

Vacuum System

A vacuum system is required to hold the thin section slides in place for sawing and grinding operations. Ward's Vacuum System (available separately) consists of a vacuum pump and a coolant trap. The black vacuum tubing should be connected as illustrated (see below)--from the pump to one flask of the coolant trap and from the second flask to the vacuum inlet at the base of the saw/grinder. The coolant trap should be checked frequently for sediment and coolant levels. If the coolant trap becomes filled with coolant and coolant is sucked into the vacuum pump, damage to the pump's components may result. To prevent damage, check coolant traps regularly and empty as needed. The vacuum may be shunted to either saw or grinder vacuum chuck by means of a valve (red ball) on the base of the unit.


**To Vacuum Intake
on Cut-Off Saw**

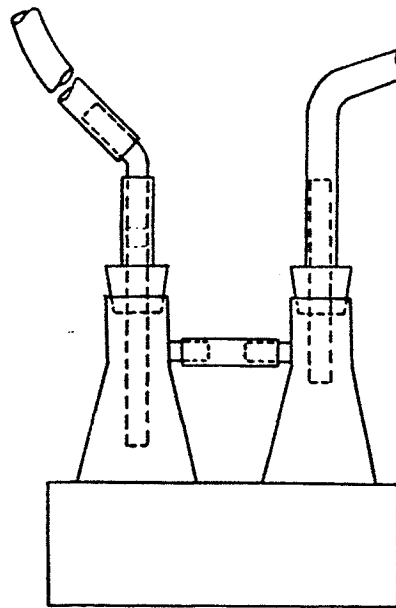

To Vacuum Pump

Coolant Trap Installation

If you plan to use both the Ingram-Ward Cut-Off Saw and separate Ingram-Ward Grinder with one common coolant and vacuum system, this may be accomplished with the set-up illustrated below. In this instance, it will be necessary to place a blank microscope slide over the openings on the vacuum chuck for the machine not in use to insure proper vacuum pressure for the machine in operation.

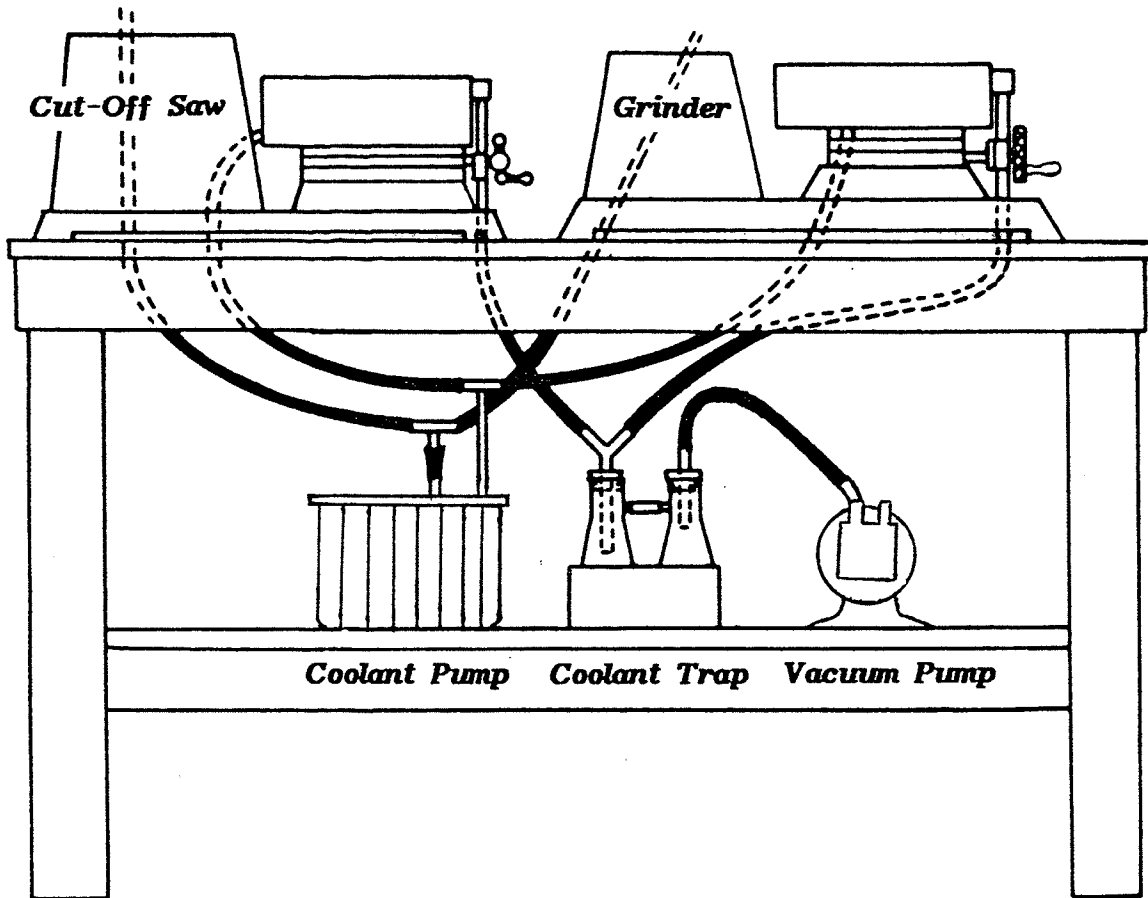

Suggested Installation Set-Up of Cut-Off Saw and Grinder using one Coolant Pump, one Coolant Trap, and one Vacuum Pump.

Electrical Hookup

Check the front of the Ingram-Ward Cut-Off Saw to make sure that all of the toggle switches on the electrical panel are in the down (off) position. At the rear of the unit, labelled electrical outlets can be found. The outlet marked VACUUM is for the vacuum pump plug and is controlled by the vacuum toggle switch on the front panel. Similarly, the outlet marked COOLANT is for the coolant pump plug and is also controlled by a single toggle switch on the front panel. These outlets should be used when a single coolant pump and a single vacuum pump system are used.

The Ingram-Ward Petrographic Thin Section Cut-Off Saw is wired for 110 volt, 60 Hz operation (unless specified for overseas shipment) and comes with a three-prong plug. It should only be plugged into a receptacle of the same voltage. **DO NOT TURN ON THIS EQUIPMENT UNTIL THE FOLLOWING CHECKLIST HAS BEEN COMPLETED.**

- (1) Check the vacuum and coolant tubing for kinks and make sure all connections are tight. Also check tubing for any sludge accumulation.
- (2) Check to see if water jets are properly adjusted with the water flinger and that they are not touching the saw blade or spindle.
- (3) Check the rocker arm assembly to make sure it has sufficient clearance from the saw blade and will not hit it if moved forward accidentally. Be sure the rocker arm is also resting on the splash shield brace before starting motor.
- (4) Be sure rubber pads are properly positioned under the corners of the machine.
- (5) Check to see that the pins on the vacuum chuck plate do not extend beyond the thickness of your microscope slide.
- (6) Empty the coolant trap in the vacuum system.
- (7) Turn on the vacuum pump.
- (8) Turn on the coolant pump.
- (9) Make sure coolant is properly directed to saw blade. Adjust pressure as needed.
- (10) Turn on the saw motor.

GENERAL OPERATING PROCEDURES

Saw Blade Preparation

It is absolutely necessary to dress (sharpen) the diamond saw blade regularly in order to keep it in proper condition for cutting. It is a simple process that requires only a few seconds but is often neglected. Please read the following instructions carefully.

Saw Blade

The cutting edge of a sintered diamond blade is made up of diamond particles in a matrix or bonding material of metal. Actual cutting is performed only along this thin rim of diamond impregnated matrix. For this surface to cut, the diamond particles must be exposed. The dressing process is one way of abrading away the metal bond with an abrasive dressing stick, furnished with the machine, thereby exposing the angular particles of diamond for better cutting. This is done quickly and easily by feeding the dressing stick into the rotating saw blade by hand and making several 1/4 inch or so cuts. Always use abundant coolant. A new diamond blade must always be dressed before attempting a cut.

For most materials, an average rate of cut should be about one square inch per minute. Softer or harder materials require less or greater time. When the rate of cutting appears to become too slow, DO NOT resort to greater pressure. This may worsen the situation. Stop cutting and re-dress the blade. Certain materials, such as very hard, compact granites, quartzites, and particularly cherts and agates, tend to burnish the metal bond over the diamond and reduce the cutting action. Forcing the blade to continue cutting under these conditions may irreversibly damage the blade. Simply dress the blade regularly when cutting seems reduced. When properly maintained, the diamond saw blade provided should last for several thousand cuts, depending upon the type of materials being sawn. If the blade ceases to cut properly even after dressing, it may be time to replace it. Inspect the rim of the blade for wear. If the diamond matrix material on the rim has been worn down to the thinner metal core of the blade, it's time for a new blade.

Making Thin Sections

Developing capability in the art of thin sectioning requires proper technique and procedure to ensure a quality end product. The key factors for making quality thin sections with the Ingram-Ward machines include proper chip preparation, uniform thickness of microscope slides, and even application of the mounting medium. When skill and technique in these areas have been achieved, the Ingram-Ward units will produce sections of excellent quality. These critical techniques of specimen preparation and mounting are explored later in this manual under "Keys to Better Thin Section Preparation". In this section, however, we will cover the general operating procedures for the cut-off saw once a specimen has been prepared and mounted for sectioning.

Cut-Off Saw Operation

The Ingram-Ward Cut-Off Saw is designed to remove the bulk of the prepared specimen chip from the slide, leaving a relatively thin layer of material for finish grinding. In the Model 65C, the saw and grinder are designed into a single unit with a common arbor, while in the Universal and Standard Ingram-Ward machines (Models 400U, 400, 137U, 135) the saw and grinder are stand alone units.

Before operating the saw:

- (1) Make sure the face plate of the vacuum chuck is completely free of dirt or grit. To clean, use a lint-free cloth.
- (2) Inspect the back surface of the specimen slide to be sawn and remove any excess mounting medium that may appear on the surface with a razor blade.
- (3) Turn on the vacuum pump. On Model 65C, make sure the vacuum switch (red ball) at the front of the unit is in the proper position for sawing. Moisten the back surface of the glass microscope slide to which the chip has been cemented. Depress the vacuum release valve and carefully place this slide surface on the face plate of the vacuum chuck. Avoid sliding your sample onto the vacuum chuck, as this may cause scratching of your slide. Make sure the microscope slide fits against all four stop pins on the chuck. At this time, it is well to check the vacuum system for possible leaks. This may be accomplished as follows: After vacuum pump has run for a short while with the microscope slide in place on the chuck, test to see if it is being held firmly in place. Leave the microscope slide in place; turn off the vacuum pump. Wait a few seconds and test again by trying to remove the slide. Now activate the vacuum release valve. If air can be heard rushing into the valve, the system is free of leaks. If, however, the slide loosens quickly when the vacuum pump is turned off, or no air rush is heard upon opening the valve, a vacuum leak is present. The most common leaks occur around the rubber stoppers in the coolant trap. Make sure they are fully seated in the flasks. Dirt or grit on the vacuum chuck or excess epoxy on the back of a slide can also cause an improper seal.
- (4) After positioning the slide on the vacuum chuck, set the micrometer on the side of the machine so that the desired thickness of the chip will be left on the slide after the saw cut. A good average thickness is .020 inches. Once the desired thickness is established, the micrometer setting should be left in this position for subsequent sawing. The graduated collar may be set at "0" for this position if desired. Remember, the glass microscope slide and the mounting medium layer will vary somewhat in thickness. This will result in a thickness variation in the sample when the micrometer is left in a fixed position. Obviously, the thicker the slide, the thinner the remaining sample after sawing.

- (5) Turn on the coolant pump and saw motor. Adjust the coolant valve so a firm spray of coolant bathes both sides of the cutting edge of the blade. It may be necessary to bend the coolant lines slightly. Direct the lines into the water flinger. In all saw models, be sure that sufficient coolant is used. Too much coolant does no harm. Too little coolant can cause excessive blade wear and possible overheating of the chip. If the blade itself overheats, some diamonds may work loose from the bonding material, thereby decreasing the life of the blade.
- (6) Fold the hinged Plexiglas guard over against the main coolant hood.
- (7) Hold the rocker arm loosely by the handle and slowly move the sample into the blade. Visually check to make sure your thickness setting leaves adequate clearance for the vacuum chuck before fully engaging. Hitting the blade with the metal stop pins or chuck will likely damage both the blade and the chuck. Once your slide sample is clearly engaged, maintain a steady uniform pressure until near the end of the cut. Toward the end of the cut, reduce the pressure slightly. The rocker arm should be moved fully forward until the stop bar is reached, then retract the rocker arm to its original position with the motor still running. When cutting, DO NOT force the sample into the blade. Apply only sufficient pressure to draw the sample at an even rate through the blade. Dress the blade as needed.
- (8) Activate the vacuum release valve and remove the slide from the chuck by lifting with your thumbnail. The sawn section is now ready to be placed on the grinder for finish grinding.

MAINTENANCE

Generally speaking, the Ingram-Ward Thin Section Cut-Off Saws are relatively maintenance-free. If kept clean and not abused, they will produce excellent thin sections for many years, requiring only periodic replacement of diamond saw blades as use demands. Follow these few basic procedures to assure quality thin section production and to help maintain equipment operating efficiency and condition:

- (1) After completing use of the equipment, wipe bare metal surfaces dry and periodically coat them with a light machine oil. DO NOT use penetrating oils, as these can be corrosive.
- (2) Periodically remove and clean the vacuum face plate pins. These can seize in place after continuous use and become difficult to remove. A pin vise is recommended for removing and replacing these pins. Replace the pins, making sure they are seated deep enough not to damage the diamond saw blade or the cup wheel during operation of the rocker arm. The pins should not extend beyond the thickness of the microscope slides in use. They should always be seated slightly below slide level so that they do not come in contact with the saw blade or cup wheel during operation.
- (3) When operating this equipment, keep coolant return hoses free of foreign material buildup. Coolant should never be allowed to accumulate in the coolant catch pan to the level of the rocker arm assembly bearings. Keep drain clear, and remove excess chip material from catch pan after use each day.
- (4) If the coolant traps overflow into the vacuum pump, simply place your finger over the exhaust nozzle of the pump, creating pressure to blow the coolant from the unit. Remove the vacuum hose from the pump during this procedure. If, after a period of time, the vacuum does not seem to be drawing air properly, some internal components may have either weakened or clogged. Repair or replacement may be required.
- (5) Keep micrometer slide advance free of dirt and grit by cleaning periodically with compressed air.
- (6) NEVER LOOSEN BOLTS on the machine (except when changing saw blades or grinding wheels), especially on the rocker arm assembly. Alignment of this assembly with the cutting surface is critical. If a suspected problem with alignment occurs, refer to the troubleshooting portion of this manual first for possible causes. If procedural steps do not correct the problem, contact Ward's Customer Service Department for technical assistance. DO NOT attempt alignment repairs without factory assistance. Most such repairs can only be done by factory-trained personnel.

TROUBLESHOOTING

Defect	Possible Cause	Remedy
Slides breaking when approaching 30 micron thickness	Different coefficients of expansion of epoxy, glass, and chip set up stresses which can cause fractures	Review mounting temperature and materials used. Equal balance.
Saw marks across chip	Chip fed into saw too rapidly	Slow feed rate
	Drifting saw blade	Dress saw blade more frequently. Reverse saw blade periodically
	Damaged saw blade	Examine blade for rim damage. Repair or replacement may be required.
Breaking slides	Micrometer feed too fast	Slow feed rate
	Saw blade drifting into glass slide	Reverse saw blade periodically.
	Loaded or damaged saw blade	Dress blade and inspect for damage.
	Vacuum not holding	Check vacuum pump intake and clean vacuum chuck plate and back of glass slide. Check hoses for leaks.
	Feeding chip into saw too fast	Slow feed rate
Chip pulls off glass slide	Micrometer feed too fast	Slow feed rate
	Improper epoxy mix (ratio of hardener to resin)	Review mixing instructions and methods
	Bonding between chip and slide inadequate	Frost the slide before mounting chip to increase surface area for greater bond.
	Air entrapped in epoxy layer	Review mounting technique
Scratches on slide	Grit, dirt, or foreign matter on vacuum chuck plate	Clean chuck plate

KEYS TO BETTER THIN SECTION PREPARATION

Introduction

Your new Ingram-Ward Thin Section equipment has been built to exacting tolerances which permit preparation of a finished thin section to a thickness of 30 microns or less without additional hand finishing in most instances. Precise, dependable equipment is obviously critical to successful thin section production. However, having the best thin section equipment does not always guarantee a quality thin section. A machine will only produce a section equal to the level of skill and technique practiced by the technician. It would be fair to say, from our experience, that 95% of the problems attributed to machine misalignment or malfunction (not grinding flat, etc.) are really the fault of improper technique. This is often the primary reason many experience inconsistent results with their thin section production. No equipment can make up for deficiencies in preparation technique. It is important for technicians to spend the necessary time at the early critical stages of preparation where their efforts will assure them a better end product.

In this section, we will review some basic procedures of thin section preparation that have proven effective in Ward's own thin section laboratory. Special emphasis will be placed on the critical areas of initial specimen grinding, microscope slide selection, and mounting.

What is a Thin Section?

In the field of Geology, a petrographic thin section is normally a fragment of rock or mineral which has been mounted on a glass microscope slide and ground to a paper thin layer (usually 0.03mm or 30 microns) for optical study. Obviously, to achieve such a thin layer requires special procedures and equipment. The majority of problems encountered during this process lie in trying to achieve a flat, thin section that is uniform in thickness and shows no evidence of major fracture or loss of specimen area. Here the technique and preparation prior to use of thin sectioning equipment is critical. All variables in microscope slide uniformity, chip or specimen flatness, and an even epoxy or other cement layer must be worked out before the final steps on the equipment. We will assume the quality of the glass slide and its index of refraction are adequate, and the mounting medium being used is a good workable product with a refractive index equal to that of the glass slide.


Specimen Preparation

The specimen to be sectioned should be washed in a suitable medium and labelled. A chip should then be cut from the bulk sample to an appropriate size for mounting using a diamond slab or trim saw. The length and width of the chip will depend upon the size of the thin section desired. For most standard petrographic slides (27 x 46 mm), a chip size of around 25 x 35 mm is usually good. The thickness of the chip should be enough to grasp easily with your fingers, but generally not more than 1/2" to 3/4" so that it will easily clear the Plexiglas coolant hood on the Ingram-Ward Cut-Off Saw when engaged. Squareness of the chip will also facilitate more efficient sawing.

If the specimen is particularly porous or exhibits many open voids, surface coating or impregnating the chip may be desired before grinding. In this case, coarse grind the surface of the chip on a flat lap and wash the chip when finished. Allow the sample to dry on a hot plate. Mix the epoxy or other mounting medium according to directions, and thin slightly with xylene or other solvent. Using a wooden spatula or other tool, apply the epoxy mixture to the ground surface of the chip until the epoxy will no longer permeate the surface. Level the epoxy by placing a glass slide over the chip and then carefully slide it off to reduce excess and eliminate grinding through an uneven surface. Cure the epoxy-coated chips on a hot plate until the epoxy is hard. If multiple sections are to be made from the sample, you may wish to impregnate the chip by immersing it in the epoxy mixture under a vacuum for 2-3 hours. Occasionally break and restore the vacuum seal to help force the impregnating solution into the specimen. Allow the chip to dry thoroughly before further grinding.

Grinding the Chip

After the chip has been cut to size and surface coated if needed, the next concern is achieving a flat chip surface for mounting. In preparing a specimen for mounting, you can use sophisticated lapping machines with diamond, or a simple glass plate with silicon carbide. Whatever you use may be dictated by cost and desired output, but the characteristics of the finished chip or specimen should be the same. After lapping with a progression of silicon carbide grits or diamond disks (to a minimum grit size of 600), the surface of the chip to be mounted must be perfectly flat and smooth. Check for surface flatness periodically with a proven technique. A "toolmaker's" knife edge can be used for this purpose. (See Diagram A.) Rest the knife edge on the ground surface of the chip while holding it up to a light source. Looking straight across the flat plane of the chip, check for light passing between the knife edge and the ground surface from length to width for any irregularities or low points. If light does appear between this contact zone in any one of the positions, the chip is not uniformly flat and should not be mounted. Lapping should continue until no variations are noted.

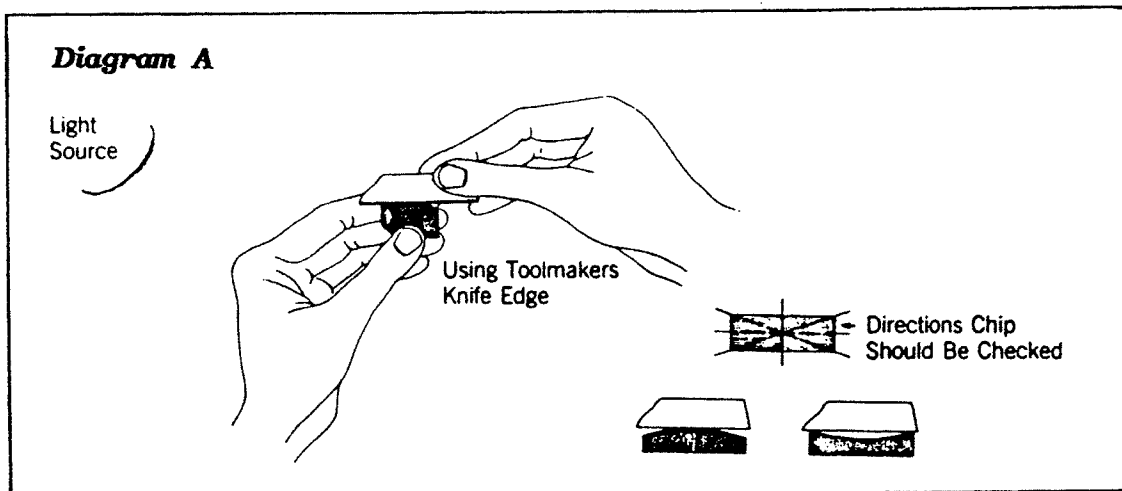

Lapping is one of the most critical areas of thin section preparation. When not performed correctly, it will result in the loss of the thin section later on during final machine grinding. (See Diagrams B & C.) A few minutes spent developing a good lapping procedure can save many hours of doing slides over again. Once this procedure has been worked out, it will become less necessary to check for flatness on each sample as you become more skilled and accustomed to your own technique and lapping ability.

Diagram B



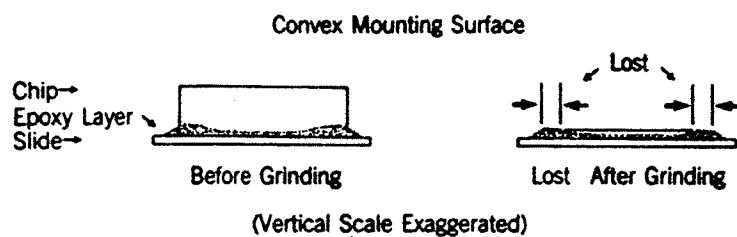


Diagram C

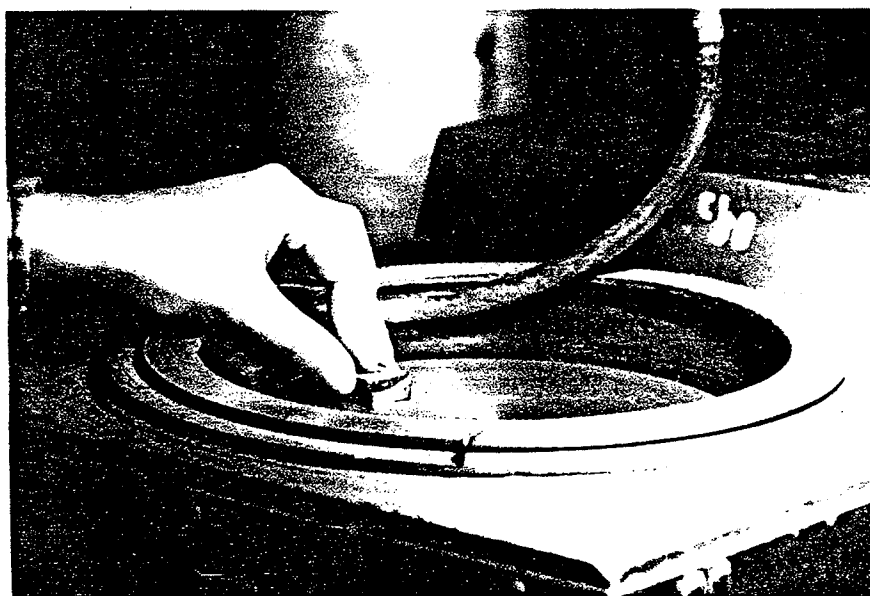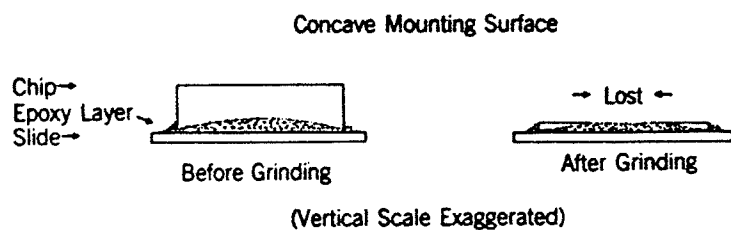

Surface Lapping of Chip

Microscope Slide Variations

The second area to consider in thin section preparation is the glass slide. Many slides will vary in thickness, from one end to the other, by more than .001 inches. This variation will make it almost impossible to get a flat, uniform thin section from any equipment used. Since the vacuum chuck on the Ingram-Ward machines and the cutting surface are parallel and have been ground to exacting tolerances, both slide surfaces must be perfectly parallel to achieve a flat finished section.

All slides should be checked with a micrometer before being used. This applies to both plain and frosted slides. Most frosted slides are either chemically frosted or sandblasted, and any irregularities in parallelism of top and bottom surfaces would remain after treatment. Using the micrometer, measure the thickness at both ends and the middle of the slide. Subtract the lowest reading from the highest reading. Any difference greater than .0005 inches will cause serious problems when grinding your section to a tolerance of .001 inches (the standard thickness of most petrographic thin sections). The usual result is a loss of specimen area during final machine grinding or the need to do excessive hand finishing to salvage the section.

To solve this problem, you could place the blank slide on the Ingram-Ward grinder and grind it down to uniform thickness before mounting the chip, or you can simply continue checking your stock of slides with the micrometer until slides of uniform thickness are found. Generally, sections on slides with slight variations of .0005 inches or less can be salvaged fairly easily by removing the slide from the grinder before fracturing or loss of specimen area occurs. The section can then be hand finished to completion by selectively grinding on a glass plate using silicon carbide #600 grit. Taking these relatively simple steps of checking slide uniformity before mounting can save hours of wasted effort and many ruined sections.

Chip Mounting

The last major area of concern is achieving an even medium (epoxy) layer between specimen and glass slide when mounting. It always seems enough of a problem mounting a chip without entrapping air bubbles in the medium. It is important to try to eliminate these bubbles, but it is only secondary to producing an even, uniform layer. The type of medium you select for this operation can save you time and effort. A number of different epoxy products, as well as Canada Balsam, are among those media in general use. Epoxies tend to form a stronger bond, are easier to use, and have a better index of refraction. Some media cure at room temperature, others require added heat, and still others cure with ultraviolet light. Your choice of medium may be dictated by desired results and the type of specimen being mounted. Working time may also be a factor. Whether hot or cold mounting the specimen, and regardless of the medium you choose, the following procedure has proven effective once the medium has been applied to the chip.

To avoid as many air bubbles as possible, position the glass slide at a 45 degree angle along the edge of the chip. (See Diagram D.) Carefully lower the glass slide until complete contact has been made between the chip surface and the slide. Work out any trapped air bubbles by applying light pressure to the back surface of the slide with the eraser end of a pencil or pointed wooden dowel. After all air is removed, press firmly in the center of the slide with the dowel (or pencil eraser) and rotate the slide in a circular motion with your other hand. (See Diagram E.) This pressure and slide movement will disperse the medium evenly and result in a uniform layer between chip and slide. (See Diagram F.) Failure to achieve this uniform layer may result in loss of specimen area during machine grinding. (See Diagram G.) Do not apply too much pressure, or the medium might be completely forced from the center, creating a dry spot. When the mount has dried thoroughly, the sample is ready for sawing and finish grinding on your thin section equipment. (See GENERAL OPERATING PROCEDURES, page 5.)

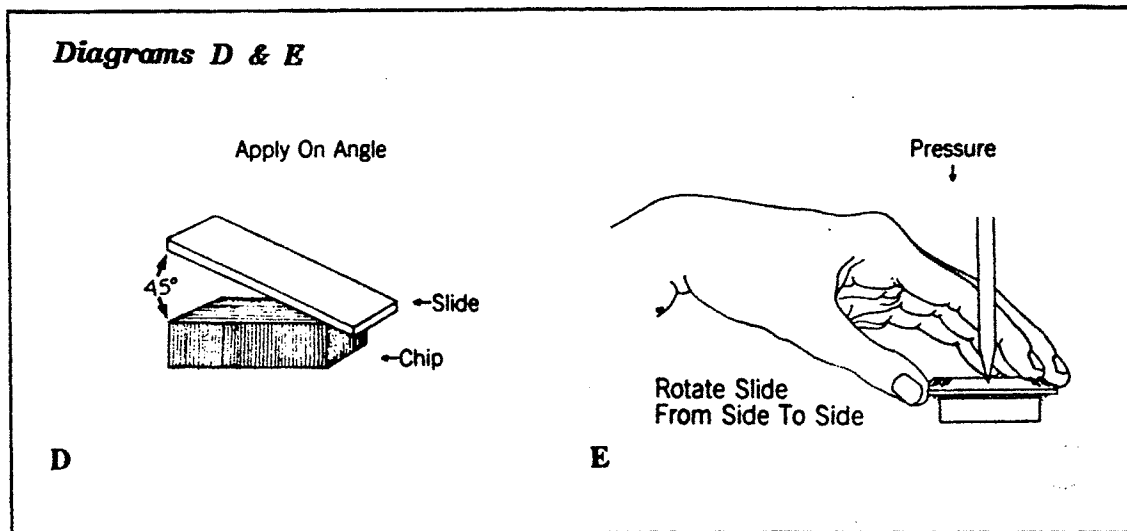


Diagram F

Flat, Smooth Chip With Even Epoxy Layer



(Vertical Scale Exaggerated)

Diagram G

Unequal Pressure



(Vertical Scale Exaggerated)

Final Preparation

The final step in completing the thin section is the application of the cover slip. The cover slip protects the finished ground surface of the sample and improves clarity for optical viewing. The cover slip should be applied with a medium having an index of refraction equal to glass and be free of impurities. After your section has been ground to final thickness, wash the section and allow it to dry in a dust-free environment. Apply a thin layer of covering medium on the section and slowly lower the cover slip onto the finished slide. Carefully remove any air bubbles as outlined under Chip Mounting. Don't apply too much pressure, or the thin cover slip will crack. After hardening, any excess media can be carefully removed with a razor blade. The section is now ready for study.

CUSTOMER SUPPORT

Your new Ingram-Ward Thin Section equipment has been designed, manufactured, and tested to perform efficiently and dependably for many years of service. The procedures and techniques outlined in this manual are designed to provide basic background information which will further insure the longevity and performance of your equipment. Please read all instructions carefully. In the event you experience difficulty in the operation of any Ingram-Ward equipment, Ward's professional staff is ready to help.

--If you are experiencing difficulty in setting up or operating the equipment, or require technical advice on proper procedure or technique, contact Ward's Thin Section Manager.

For Technical Assistance CALL TOLL FREE: 1-800-962-2660.

--In the event repairs or replacement parts are ever needed, all work is handled through our manufacturer in Griffin, Georgia. Contact Ward's Customer Service Department for details and shipping instructions.

For Customer Service CALL TOLL FREE: 1-800-962-2660.

SUPPORT EQUIPMENT

The following support equipment is available separately from Ward's and is designed for use with all Ingram-Ward Thin Section equipment. All Ingram-Ward saws and grinders require a coolant system, coolant trap, and vacuum pump for operation. Suitable equipment may already be available in your laboratory, or you may wish to purchase these specific Ward's products. Complete descriptions and pricing information may be found for these and other related products in your Ward's Geology/Earth Science Catalog.

Catalog Number	Description
28 E 1920	Coolant Pump System
28 E 1921	Coolant Trap
21 E 7410	Vacuum Pump
15 E 8570	Mobile Thin Section Lab Table
28 E 1901	Replacement Diamond Saw Blade - 5 inch (For Models 65C and 135)
28 E 1907	Replacement Diamond Saw Blade - 7 inch (For Model 137-U)
28 E 1915	Replacement Diamond Cup (Grinding) Wheel (For all Models 65C, 400, and 400-U)
28 E 1927	Dressing Sticks for Cut-Off Saws (1" x 1" x 6")
28 E 1928	Dressing Sticks for Grinders (1/2" x 1/2" x 6")

Accessory Materials - Suggested items for basic thin section production needs.

Catalog Number	Description
25 E 4600	Polarizing Film Disks - 2-inch diameter, set of 2. (For checking specimen thickness/birefringence.)
15 E 4700	Metric Micrometer (For checking slide/specimen thickness.)
13 E 0801	Petrographic Microscope Slides, Etched 27 x 46 mm, Package of 72
13 E 0800	Petrographic Microscope Slides, Plain 27 x 46 mm, Package of 72
14 E 3245	Cover Slips (glass) - 24 x 40 mm, 1 oz. package

Accessory Materials - continued

Catalog Number	Description
15 E 4509	0151 Epoxy Mounting Medium - Kit, 2 tubes (Resin - 2.54 oz./Hardener - .81 oz.) Sufficient for mounting several hundred slides.
37 E 9550	CR-1 Cover Glass Mounting Medium - 4 oz. bottle.
	Silicon Carbide Abrasives
28 E 1480	80 Grit
28 E 1490	100 Grit
28 E 2510	120 Grit
28 E 2520	220 Grit
28 E 2530	320 Grit
28 E 2540	400 Grit
28 E 2560	600 Grit
14 E 3582	Slide Warmer Table - 0 to 100 degrees C. (For hot specimen mounting and curing.)
28 E 1233	Petrographic Specimen Polisher - 5 speeds, 115V. (For lapping chips.)
28 E 1658	Slab and Trim Saw (10") - 110V with diamond blade. (For cutting and sizing chips.)
30 E 4801	Petrographic Slide Storage Box Capacity: 100 slides, 27 x 46 mm.

WARD'S

Ward's Natural Science Establishment, Inc.
5100 West Henrietta Road • P.O. Box 92912
Rochester, New York 14692-9012 • (716) 359-2502